

ProteoSpin™ Endotoxin Removal Micro Kit
Product # 22800

Product Insert

The ProteoSpin™ Endotoxin Removal Micro Kit is designed for the rapid removal of endotoxins from up to 150 µg of previously purified proteins or peptides. Endotoxins, also known as lipopolysaccharides, are cell-membrane components of Gram-negative bacteria such as *E. coli*. Endotoxins liberated by Gram-negative bacteria are frequent contaminations of protein solutions derived from bioprocesses. Due to the high toxicity of endotoxins *in vivo* and *in vitro*, their removal from protein preparations is often necessary prior to the use of the protein in downstream applications. The ProteoSpin™ Endotoxin Removal Micro Kit efficiently reduces endotoxin levels to ≤ 0.01 EU/µg of protein, using spin column chromatography based on Norgen's proprietary resin as the separation matrix. The purified protein samples can be used in a number of downstream applications including sequencing, cloning, and *in vitro* and *in vivo* introduction into cells and organisms for various purposes.

The ProteoSpin™ Endotoxin Removal Micro Kit contains sufficient materials for 20 purifications. Each column is able to remove endotoxins from up to 150 µg of previously purified proteins, with protein recoveries of > 95% being achieved. Preparation time for a single sample is approximately 20 minutes.

Specifications

Kit Specifications	
Maximum Protein Input	150 µg
Maximum Protein Volume Input	450 µL
Final Endotoxin Levels	≤ 0.01 EU/µg protein
Average Recovery	>95 %
Minimum Elution Volume	50 µL
Time to Complete 10 Purifications	20 minutes

Kit Components

Component	Product # 22800 (20 samples)
Binding Buffer	4 mL
Activation and Wash Buffer	50 mL
Elution Buffer	6 mL
Endotoxin Removal Solution	1.5 mL
Neutralizer	2 mL
Spin Columns	20
Collection Tubes	20
Elution Tubes (1.7 mL)	20
Product Insert	1

Storage Conditions and Product Stability

All solutions should be kept tightly sealed and stored at room temperature. All the reagents should remain stable for at least 1 year in their unopened containers.

Precautions and Disclaimers

This kit is designed for research purposes only. It is not intended for human or diagnostic use.

Ensure that a suitable lab coat, disposable gloves and protective goggles are worn when working with chemicals. For more information, please consult the appropriate Material Safety Data Sheets (MSDSs). These are available as convenient PDF files online at www.norgenbiotek.com.

Customer-Supplied Reagents and Equipment

- Benchtop microcentrifuge
- Microcentrifuge tubes
- Isopropanol
- Sterile, deionized water or Milli-Q[®] water

Procedure

All centrifugation steps are carried out at 14,000 x g in a benchtop microcentrifuge. Please check your microcentrifuge specifications to ensure proper speed. Performance of the kit is not affected by temperature, and thus the procedure may be performed at room temperature, 4 °C, or on ice.

Notes prior to use:

- Ensure that all solutions are at room temperature prior to use, and that no precipitates have formed. If necessary, warm the solutions and mix well until the solutions become clear again.
- Ensure that the maximum protein input does not exceed 150 µg or 450 µL. If the mass or volume of protein exceeds this, the sample will need to be processed using more than 1 column.
- High molarities of salts will interfere with the pH adjustment in the Sample Preparation step. Dilute the molarity of the input protein solution to < 50 mM using sterile, deionized water or Milli-Q[®] water. If the volume increases over the maximum 450µL input, the sample can be split and processed using 2 columns.

1. Column Activation

- a. Open the cap on the pre-assembled spin column with its 2 mL collection tube.
- b. Apply 500 µL of **Column Activation and Wash Buffer** to the column and close the cap.
- c. Centrifuge for 1 minute and discard the flowthrough.
- d. Repeat steps **1b** and **1c** to complete the column activation step.

2. Sample Preparation

- a. Obtain protein sample and measure the volume. Bring the volume up to 450 µL using sterile, deionized water or Milli-Q[®] water.
- b. Add 20 µL of **Binding Buffer** to the protein sample and mix well.
- c. Verify that the pH is 3.5 or below, and add more pH Binding Buffer if necessary in 20 µL increments.
- d. Transfer the protein sample into the top reservoir of the activated column (from Step 1 above).
- e. Add a 1% volume of **Endotoxin Removal Solution** to the liquid on top of the column. (i.e.: Add 5 µL for every 500 µL volume of input placed onto the column).

Note: The Endotoxin Removal Solution is extremely viscous, and thus care should be taken when measuring out the correct volume.

- f. Close the lid of the column and vortex the column assembly **GENTLY** to mix. Let stand for 5 minutes at room temperature.

Note: Some dripping of the protein sample into the collection tube may occur during the incubation period. Even if this does occur, proceed with the protocol as written.

- g. After the 5 minute incubation, add a 10% volume of **isopropanol** (supplied by the user) to the liquid on the top of the column. Close the lid and vortex column assembly **GENTLY** to mix. (i.e.: Add 50 μL of isopropanol for every 500 μL of sample on top of the column).

2. Binding to Column

- a. Place the column into a microcentrifuge and spin at 14,000 x g for 1 minute.
- b. Discard the flowthrough and reassemble the spin column with its collection tube.

3. Washing Bound Protein

- a. Apply 500 μL of **Activation and Wash Buffer** to the column assembly and centrifuge the unit for 1 minute at 14,000 x g.
- b. Discard the flowthrough and reassemble the spin column with the collection tube.
- c. Repeat steps **3a** and **3b**.
- d. Spin the column for an additional 1 minute at 14,000 x g, in order to completely dry the resin. Discard the collection tube.

4. Elution of Clean Protein

- a. Add 9.3 μL of Neutralizer to a provided **Elution tube**.
- b. Assemble the column (with protein bound to the resin) with the **Elution tube** from above.
- c. Add 50 μL of **Elution Buffer** to the center of the resin bed and centrifuge the column assembly for 1 minute at 14000 x g.
- d. Without removing the assembly, add another 50 μL of **Elution Buffer** and centrifuge for 1 minute at 14,000 x g to elute the proteins.

Note: The majority of the input proteins will be recovered in the first elution. However, a second elution may be performed if desired. Steps 4a to 4c should be repeated, and the elution should be collected into a fresh elution tube in order to prevent dilution of the first elution. The total elution volume can also be reduced to 50 μL if required (1 x 50 μL elution), but the amount of Neutralizer must be reduced to 4.7 μL as well.

Troubleshooting Guide

Problem	Possible Cause	Solution and Explanation
Protein solution does not flow through the column	Centrifugation speed was too low	Check the centrifuge and ensure that it is capable of generating 14,000 x g. Sufficient centrifugal force is required to push the liquid through the column.
	Inadequate spin time	Spin for an additional minute to ensure that the liquid is able to flow through the column.
	Cellular debris is present in the protein solution	Prior to the sample preparation step, filter the sample with a 0.45 µM filter or spin down insoluble materials. Solid, insoluble materials can cause clogging problems.
Poor Protein Recovery	Input protein solution is very high in molarity.	The Binding Solution may not lower the pH of the protein solution sufficiently if the input protein solution has a very high molarity. Dilute the molarity of the input to < 50mM with water. If the volume of the input amount is increased over the maximum 450 µL input, the sample can be split and processed using 2 columns.
	The appropriate amount of Binding Buffer was not added	Ensure that 20µL of Binding Buffer is added for every 500 µL of protein processed. The protein volume must not exceed 450 µL.
Protein does not perform well in downstream applications	A different Elution Buffer was used	The provided Elution Buffer has been optimized for endotoxin-free recoveries. The endotoxin-free properties of the eluted protein will be compromised if another elution buffer is used. If a different Elution buffer other than the one provided is used, the buffer should also be checked for any components that may interfere with the application. Common components that are known to interfere are high salts, detergents and other denaturants. Check the compatibility of your elution buffer with the intended use.
Endotoxin levels are high in the elution	A different Elution Buffer was used	The provided Elution Buffer has been optimized for endotoxin-free recoveries. The endotoxin-free properties of the eluted protein will be compromised if another elution buffer is used. If a different Elution buffer other than the one provided is used, the buffer should also be checked for Endotoxin levels.
	Extremely high endotoxin levels in input protein	Check the endotoxin levels of your protein input. The endotoxin levels of the eluted protein may be reduced significantly compared to the input.

Problem	Possible Cause	Solution and Explanation
Eluted protein is degraded	Eluted protein was not neutralized	Add 9.3 μ L of Neutralizer for every 100 μ L of eluted protein in order to adjust the pH to neutral. Some proteins are sensitive to high pH, such as the elution buffer.
	Bacterial contamination of protein solution	The protein samples can be prepared with 0.015% sodium azide.
	Eluted protein was not neutralized quickly enough	If the eluted protein is not neutralized immediately, degradation will occur. We strongly recommend adding Neutralizer in order to lower the pH.

Related Products	Product #
ProteoSpin™ Endotoxin Removal Maxi Kit	22200
Endotoxin Removal Kit (Mini)	21800
Endotoxin Removal Kit (Maxi)	21900
Plasmid MaxiPrep Kit (Endotoxin Free)	15300
ProteoLadder 100, Lyophilized	13500
ProteoLadder 100	12300

Technical Support

Contact our Technical Support Team between the hours of 8:30 and 5:30 (Eastern Standard Time) at (905) 227-8848 or Toll Free at 1-866-667-4362.

Technical support can also be obtained from our website (www.norgenbiotek.com) or through email at techsupport@norgenbiotek.com.

3430 Schmon Parkway, Thorold, ON Canada L2V 4Y6
 Phone: (905) 227-8848
 Fax: (905) 227-1061
 Toll Free in North America: 1-866-667-4362